## Colcemid Solution

Catalog No. 9311

10 mL

10 µg/mL in Hank's Balanced Salt Solution  
10 µg/ml in Hank's Ausgeglichenen Salzlösung  
10 µg/ml in Soluzione di Hank di Sale Equilibrato  
10 µg/ml en Solución Salina Equilibrada de Hank's  
10 µg/ml dans la Solution Sa line Équilibrée de Hank  
10 µg/ml em Solução Salina Equilibrada Hank's  
10 µg/mL σε διάλυμα αλάτων Hank (Hanks' Balanced Salt Solution/HBSS)  
10 µg/ml v Hankově vyváženém solném roztoku  
10 µg/ml i Hank's afbalancerede saltopøsning  
10 µg/ml Hank's Balanced Salt Solution -suolaliuoksessa  
10 µg/ml Henka līdzsvarotā sāls šķīdumā  
10 µg/ml in Hanks gebalancerde zoutoplossing  
10 µg/ml w zbalansowanym roztworze soli Hanka  
10 µg/ml in Soluție Salină Echilibrată Hank  
10 µg/ml i Hanks balanserade saltlösning  
10 µg/ml Hanki tasakaalustatud soolalahuses  
10 µg/ml Hank-féle kiegyenlített sóoldatban  
Henko subalansuotas druskų tirpalas, 10 µg/ml  
Hanks Dengeli Tuz Solüsyonunda 10µg/ml  
10 µg/ml v Hankovom vyváženom solnom roztoku  
10 µg/ml в балансиран солев разтвор на Ханк  
10 µg/ml u Hankovoj izbalansiranoj otopini  
10 µg/mL f'Soluzzjoni Salina Bilanċjata ta' Hank  
10 µg/ml v Hankovi uravnoteženi solni raztopini

## Glossary of Symbols\*:

Catalog Number

Lot Number

Sterilized using aseptic processing techniques (filtration)

Expiration:  
Year - Month - Day

Caution, consult accompanying documents

Consult instructions for use

Storage Temperature  
2-8°C

Manufacturer

CE Mark

Emergo Europe - Westervoortsedijk 60  
6827 AT Arnhem  
The Netherlands\*Symbol Reference - EN ISO 15223-1, Medical devices –  
Symbols to be used with medical device labels, labeling.

## REFERENCES

1. L.S. Goodman and A. Gilman. *The Pharmacological Basis of Therapeutics*. 5<sup>th</sup> edition. p. 351. (Mac Millan. New York. 1975).
2. T. T. Puck. *Studies of the Life Cycle of Mammalian Cells*. Cold Spring Harbor Symposia. Vol. 29. pp. 167-176. (1964).
3. J. J. Yunis. *Human Chromosome Methodology*. 2<sup>nd</sup> edition. p. 186. (Academic Press. New York. 1974).
4. M. H. Wolff and K. E. Schneewies. *The Multiplication of Herpes Simplex Virus in the Different Phases of the Cell Cycle of Hep-2 Cells*. Zbl. Bakt. Hyg., I. Abt. Org. A. 224. pp.145-155.
5. E. Robbins. *Cell Synchronization in Methods in Enzymology*. S. P. Colowick, N. O. Kaplan Ed. Vol XXXII Biomembranes Part B. p. 596. (1974).
6. W. Nultsch. *Allgemeine Botanik* Goerg Thieme Verlag. Stuttgart, 6<sup>th</sup> edition. p. 328. (1977).
7. H. E. Street. *Single Cell Clones in Street H*. Plant Tissue and Cell Culture. P. 203. (Blackwell Oxford. 1973).

FUJIFILM Irvine Scientific, Inc.  
2511 Daimler Street, Santa Ana, California 92705 USA  
Telephone: 1 949 261 7800 • 1 800 437 5706 Fax: 1 949 261 6522 • www.irvinesci.com

© 2023 FUJIFILM Irvine Scientific, Inc. All rights reserved. FUJIFILM Irvine Scientific, and the FUJIFILM Irvine Scientific logo, are trademarks of FUJIFILM Irvine Scientific, Inc. in various jurisdictions. All other trademarks are the property of their respective owners

PN 40129 Rev. 11

Effective Date: 27-DEC-2023

## INDICATION FOR USE

Colcemid Solution is intended for *in vitro* diagnostic use.

## DEVICE DESCRIPTION

Colcemid Solution, is a ready-to-use solution with a concentration of 10 µg/mL of N-deacetyl-N-methyl colchicine (C<sub>21</sub>H<sub>25</sub>NO<sub>3</sub>) in Hank's Balanced Salt Solution. Colcemid is used to arrest mitosis in plant and animal cells by preventing spindle formation during metaphase of the cell cycle.<sup>1</sup> This effectively allows for the optimal yield of metaphases through enabling synchronization of cell cycles to facilitate cytogenetic analysis. Colcemid Solution is at least 30-40 times less toxic than colchicine.<sup>2</sup>

## QUALITY ASSURANCE

A Safety Data Sheet and lot-specific Certificate of Analysis are available upon request.

## SUGGESTED USES

1. Chromosome Analysis – amniotic cells after initial cultivation.<sup>3</sup>  
Colcemid Solution arrests mitotic cultured cells in metaphase. Karyotyping for chromosome analysis is greatly facilitated when large numbers of metaphase cells are present in culture.
2. Cell Synchronization  
Colcemid Solution aids in determining the mitotic activity index for cell synchronization studies.<sup>4</sup> Through investigations, Colcemid Solution has been shown to inhibit HeLa and CHO cells at concentrations of 0.05 – 0.5 µg/mL.<sup>5</sup>
3. Polyploid Plants and Asexual Reproduction  
Colcemid Solution produces an increase in the chromosome count. New plant species can be produced in this manner under suitable conditions.<sup>6</sup> In asexual plant reproduction, the chromosome complement can be doubled under certain conditions.<sup>7</sup>

## STORAGE INSTRUCTIONS AND STABILITY

Store tightly closed container at 2 – 8° C, protected from light.

When stored as directed, Colcemid is stable until the expiration date shown on the bottle label.

## PRECAUTIONS AND WARNINGS

Colcemid Solution is mutagenic, tumorigenic, embryotoxic and teratogenic.

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